


Ellipsometry is a powerful analytical tool in the characterisation of thin films in many applications, including semiconductors, dielectrics, metals and polymers. It is a non-contact, non-destructive optical technique, which measures the polarization change of reflected light after interaction with a layer. This change in polarisation is related to thin film properties like thickness and refractive index.

# Spectroscopic Ellipsometry

- layer thickness
- refractive index
- extinction coefficient
- multi-layers
- temperature dependence

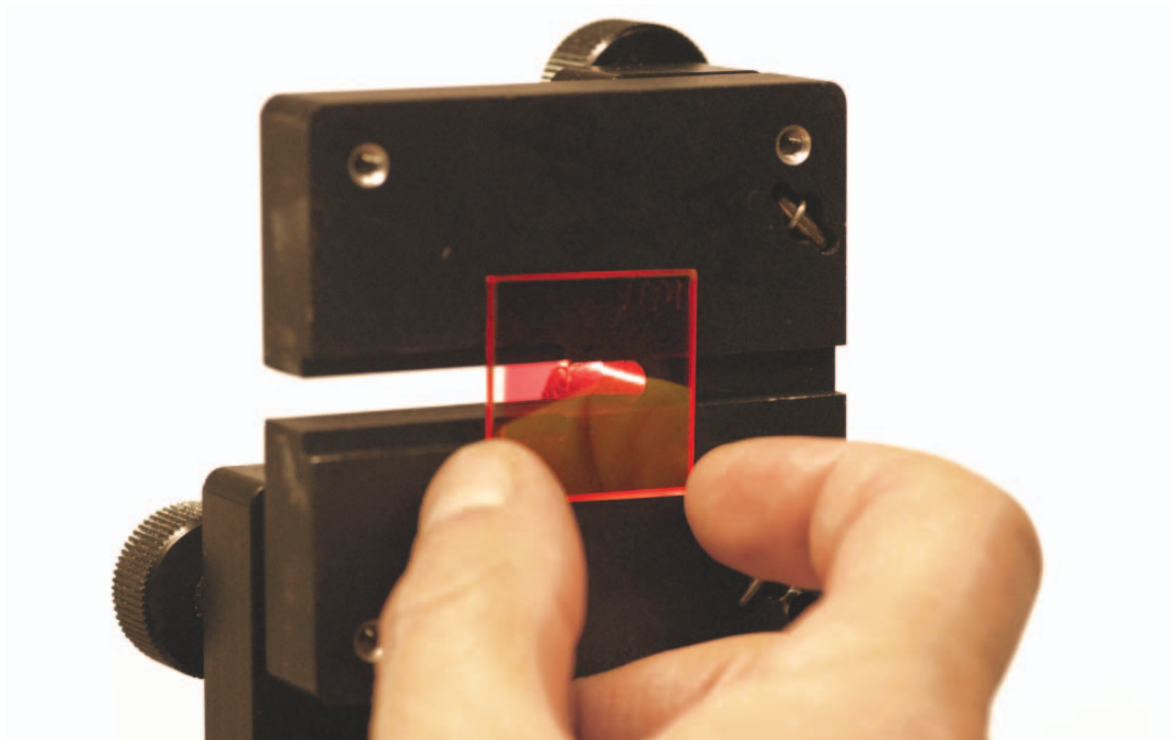

MiPlaza



**PHILIPS**

**Theory**

Ellipsometry involves light reflecting from a surface of interest. The plane of incidence is the plane perpendicular to the surface of the sample (see Figure 1). The polarization state of the light is defined by two components, one parallel (p) to the plane of incidence and one perpendicular (s).

It is important to understand that these components reflect differently from the sample, depending on the angle of incidence. Notice the differences in reflection of the "p" and "s" component in the silicon example of

Figure 2a. The  $R_s$  reflection increases steadily, while the  $R_p$  reflection goes through a minimum. This angle is called the Brewster angle. Ellipsometry measurements are most sensitive to film characteristics at this Brewster angle. As a light wave is propagating it appears to be a sinusoid from the side (Figure 3a). However, if you were to look down the z-axis with the wave coming directly into your eye, the electric field of the wave would appear to be a vertical line along the y axis as shown in Figure 3b. The polarization of a wave can be rotated to any angle. To describe its polarization, the wave

can be projected onto the X and Y axes. When a wave is defined like this, there are two electric field components,  $E_x$  and  $E_y$ . When these components are propagating in the same direction, perpendicular and in phase with each other, a linearly polarized wave results. Elliptically polarized light is illustrated in figure 4. In this polarization state, the vectors  $E_x$  (the s component) and  $E_y$  (the p component) can take on any arbitrary phase and amplitude. The resulting concept is an ellipse propagating in space.

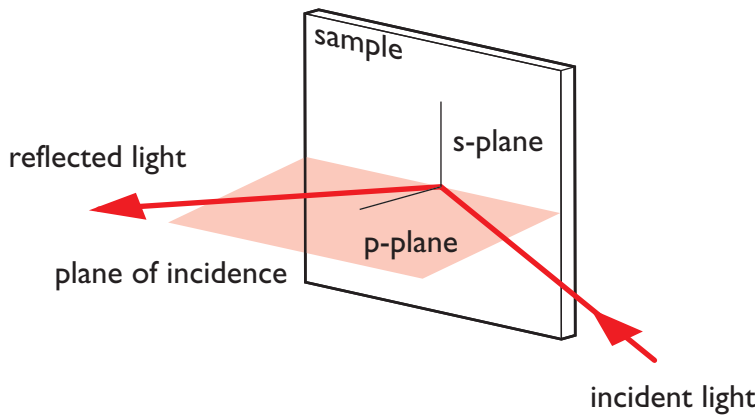



Fig 1: The plane of incidence.

**Ellipsometry**

The term ellipsometry comes from measuring elliptically polarized light. Figure 1 illustrates the basic principle behind ellipsometry. First, the polarization state of incoming light is known. This incident light interacts with the sample and reflects from it. The interaction of the light with the sample causes a polarization change in the light, from linear to elliptical polarization. This change is measured by analysing the light reflected from the sample.

To describe polarization there are two parameters, orientation and ellipticity. They are related to the two ellipsometric values, Psi ( $\Psi$ ) and Delta ( $\Delta$ ), which describe the polarization state of reflected light after interaction with a film (See Figure 2b). Because we measure a ratio and a phase change of one and the same light beam, both parameters can be measured very accurately. Measurements are not affected by light source intensity fluctuations.

Fig 2: a) the reflection as a function of angle of incidence for the parallel ( $R_p$ ) and perpendicular ( $R_s$ ) components of polarised light . b) behaviour of Psi ( $\Psi$ ) and Delta ( $\Delta$ ), which describe the polarization state of reflected light after interaction with a film.  $\Psi$  is the Ratio of the reflected intensities  $R_p$  and  $R_s$  and  $\Delta$  is the relative phase change of these two waves.

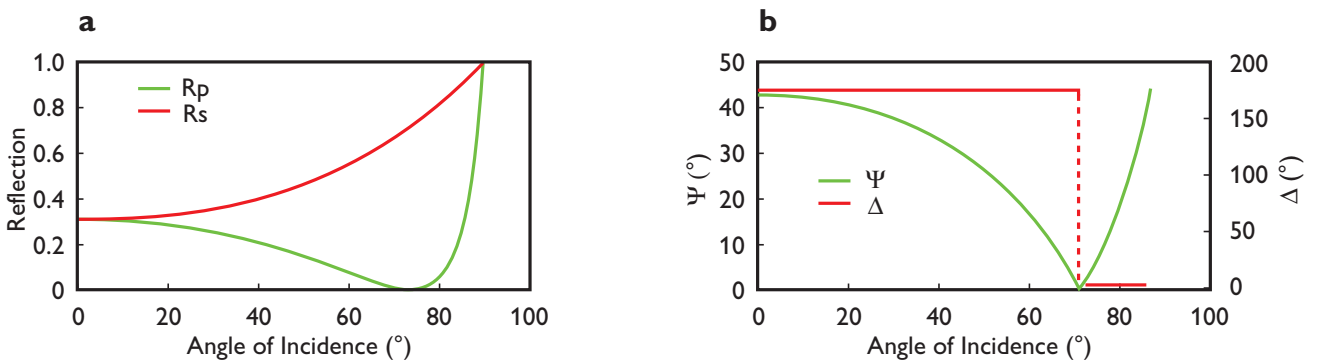

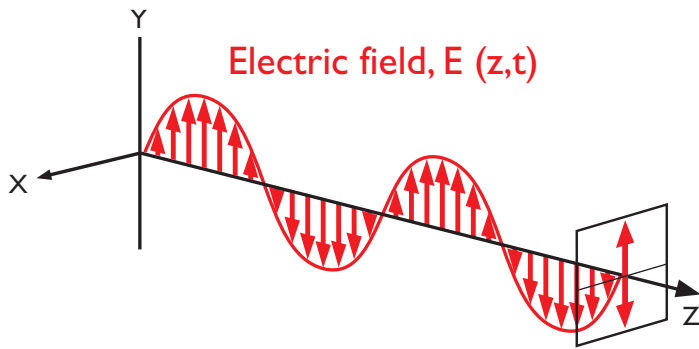


Fig. 3: Visualisation of propagating linearly polarised light.

**VASE**

Variable Angle Spectroscopic Ellipsometry (VASE) is used to measure at different angles of incidence and in a wavelength range of interest. This helps to ensure optimum sensitivity to any problem and determine multiple unknown parameters in difficult, multi-layer samples.

**Refractive index and extinction coefficient**

Ellipsometry is primarily used to measure film thickness, the refractive index (n) and the extinction coefficient (k). Both n and k are needed to describe real materials and are not constant, but vary with wavelength and temperature.

**Layer thickness**

When light is incident on a thin film it is reflected by both the top and bottom interfaces. Each reflected wave will have its own phase and amplitude, which causes interference in the reflected light. Thicker films will have more oscillations versus wavelength (Figure 5). This enables very accurate layer thickness measurements, both for single layer thin films (see Figure 6) as well as for multilayer stacks.

**Data Analysis**

Ellipsometry does not actually measure optical constants or film thickness, but Psi and Delta are functions of these characteristics. To extract useful information about thin films, a model is constructed that describes the optical parameters of the sample. Then, the unknown parameters and the thickness are fit to obtain a best match between the theoretical response and the experimental data.

Fig. 4: Visualisation of propagating elliptically polarised light. For an elliptically polarized wave both components have arbitrary amplitude and phase.


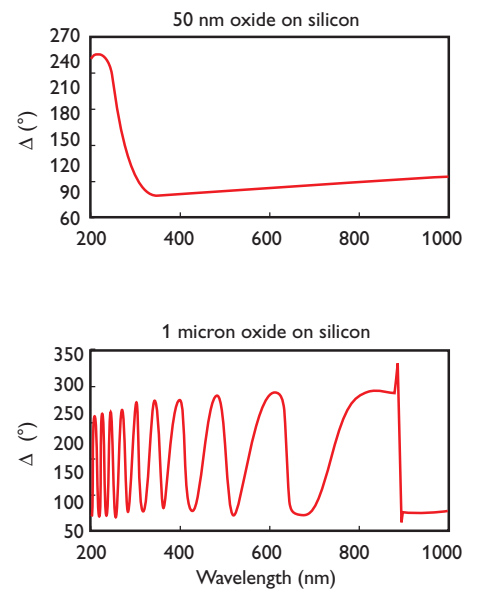


Fig. 5: Layer thickness studies: the number of oscillations is characteristic for the film thickness.

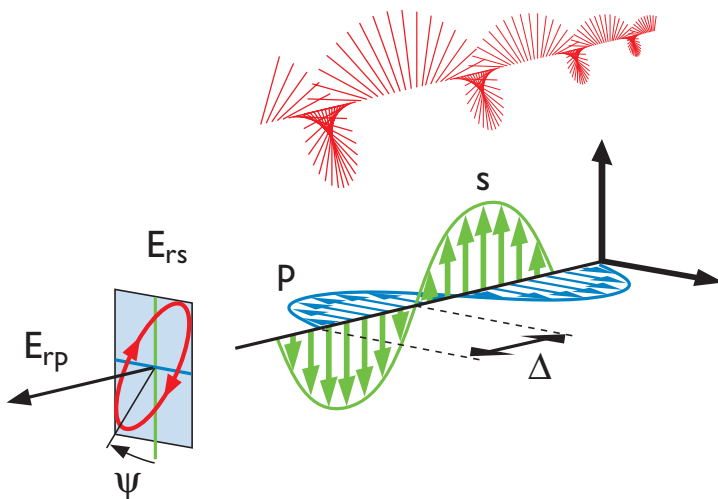

## Philips Research Materials Analysis

offers a full range of analytical methods and expertise to support both research and manufacturing, serving customers by taking an integral, solution-oriented approach.

## making the invisible visible

### For more information:

Tel./fax: +31 -40-27 48044/43075

E-mail: [materialsanalysis@philips.com](mailto:materialsanalysis@philips.com)

<http://www.miplaza.com/materials.html>

<http://pww.natlab.research.philips.com:25222/>

### Technical Note 21

May 2007

Toon de Win

### Applications

- thickness, optical constants of organic and inorganic thin films on flat substrates
- individual layer thicknesses of complex multi-stack samples
- thermal expansion of organic and inorganic thin layers
- optical anisotropy of polymer films

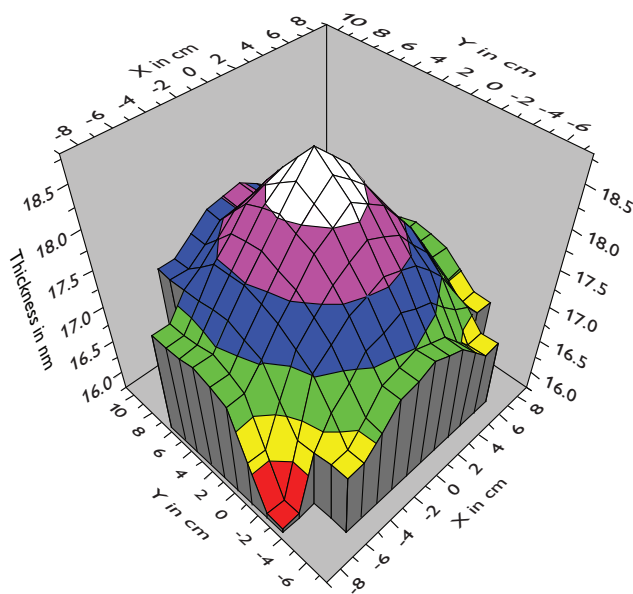


Fig. 6: Example of thickness mapping of a layer of phase change material as used as the data storage layer in DVD+RW discs.

### Characteristics

#### Measurement

Measurement of Psi, Delta data at different angles

of incidence over a wide wavelength range (193 – 1700 nm) with a variable step-size.

#### Acquired data

Reflection and transmission ellipsometric data, polarized transmission and reflection intensity, anisotropic measurements.

#### Lateral information

>1 mm, depending on spot size.

Automated sample translation stage (150 x 150 mm) available for mapping procedures.

#### Accuracy (thickness, refractive index)

Strongly dependent on sample geometry.

#### Sample type & requirements

Solids, thin layers and multilayers with optically flat substrates. Thickness from sub-monolayer coverage to  $\mu\text{m}$  range.

#### Heating experiments

Temperature (max. 300°C) dependent measurements in different atmospheres, used for the determination of the thermal expansion of thin films.

#### Spot size

200  $\mu\text{m}$  or 2 mm.

